visocolor® alpha

# Ammonium Ammonio / Amonio 0.2–3 mg/L NH<sub>4</sub><sup>+</sup>

REF 935 012

50 Tests

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## Test kit for performing colorimetric tests on ammonium

en

**Hazard warning:** NH<sub>4</sub>-1 contains sodium hydroxide solution 2–5%.

H314 Causes severe skin burns and eye damage.

**Instructions for use:** *also refer to the pictogram on the inside*

1. Rinse the test vessel several times with the water sample and fill to the ring mark (5 mL).
2. Add 10 drops of NH<sub>4</sub>-1 and mix by swirling.
3. Add 1 level measuring spoon of NH<sub>4</sub>-2 and dissolve by swirling. Wait for 5 min.
4. Add 4 drops of NH<sub>4</sub>-3 and mix by swirling.
5. After 5 min place the measuring vessel on the color chart and assign the value by comparison of the color. Mid-values can be estimated.

This method can be also used for analyzing sea water after dilution (1+9).

**Disposing of the sample:** The used analysis specimens can be flushed down the drain with tap water and channelled off to the local sewage treatment works.

**Interferences:** Strongly acidic or buffered solutions have to be adjusted to pH 9–10 for the test using sodium hydroxide solution. Primary amines react in the same way as ammonium ions and produce higher results. Depending on their concentration, substances which draw on the chlorine may reduce the measurement reading or suppress the reaction totally.

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**en** P260, P280, P301+330+331, P303+361+353, P304+340, P305+351+338, P501 Do not breathe vapors. Wear protective gloves/eye protection. IF SWALLOWED: Rinse mouth. Do NOT induce vomiting. IF ON SKIN (or hair): Take off immediately all contaminated clothing. Rinse skin with water/shower. IF INHALED: Remove person to fresh air and keep comfortable for breathing. IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing. Dispose of contents /container to regulated waste treatment. For further information ask for a safety data sheet.

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MACHERY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6–8 · 52355 Düren · Germany